

WiFi - receive side.
- glass cabinet.
- Ethernet plug

- red ethernet
cable to bottom
right port.

LASER INTERFEROMETER GRAVITATIONAL WAVE OBSERVATORY
- LIGO -

CALIFORNIA INSTITUTE OF TECHNOLOGY
MASSACHUSETTS INSTITUTE OF TECHNOLOGY

CR: 202.

LIGHTS: - Turn off in Control room.

Technical Note	LIGO-T1500062-v8	2019/06/04
Pcal End Station Power Sensors Responsivity Ratio Measurements Procedures and Log		
S. Banagiri, N. Lecoecuche, S. Karki, R. Savage		

Upon arrival: - call 202 and say we are
- key on light transitioning EX to hazard & change
in bathroom. door sign.

California Institute of Technology
LIGO Project, MS 18-34
Pasadena, CA 91125
Phone (626) 395-2129
Fax (626) 304-9834
E-mail: info@ligo.caltech.edu

Massachusetts Institute of Technology
LIGO Project, Room NW22-295
Cambridge, MA 02139
Phone (617) 253-4824
Fax (617) 253-7014
E-mail: info@ligo.mit.edu

LIGO Hanford Observatory
Route 10, Mile Marker 2
Richland, WA 99352
Phone (509) 372-8106
Fax (509) 372-8137
E-mail: info@ligo.caltech.edu

LIGO Livingston Observatory
19100 LIGO Lane
Livingston, LA 70754
Phone (225) 686-3100
Fax (225) 686-7189
E-mail: info@ligo.caltech.edu

END STATION: DATE:

PEOPLE DOING THE WORK:

Before (or after) going to the End Station

- Check the calibration of the Keithley Model 2100 voltmeter using the Martel Calibrated Voltage source (the same one that will be taken to the end station). Note: use negative polarity.
 - 4 V with Martel = V on Keithley 2100 DVM
 - 2 V with Martel = V on Keithley 2100 DVM
 - 0 V with Martel = V on Keithley 2100 DVM


Items to take to the end station for the measurements:

- Working Standard
- PD Satellite Box (blue box), D1300368
- Long (25') and short (6'-10') 9-pin D-sub cables
- BNC cable
- Martel calibrated voltage source, banana-to-BNC adapter cable, and charger/AC adapter
- IR-only laser glasses (for use ONLY if work in the green, ALS laser enclosure will not be ongoing in parallel)

Before starting the measurements

- Disable excitations on the Pcal MEDM screen. (sitemap - CAL)
- Check that ETM is pointed properly. Check Pcal beam locations at the Rx sensor (photograph spot locations on white card). X-arm ETM - ETMX
- Open gpscloc - ALIGNED TOP SCREEN
- Open StripTool and display the following sensor outputs. Always verify that signals are stable before recording time series.
 - (IFO):CAL-PCAL(END)-TX_PD_VOLTS_OUTMON
 - (IFO):CAL-PCAL(END)-RX_PD_VOLTS_OUTMON
 - (IFO):CAL-PCAL(END)-WS_PD_OUTMON
 - (IFO):CAL-PCAL(END)-OFS_PD_OUTMON

} dog from sitemap.

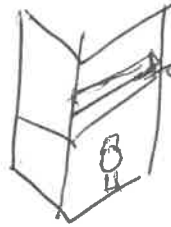

short DB9
from power box
back
(NO BNC
on back).

Calibration measurements and log

Preliminary measurements

Calibrate the Working standard channel


- Connect Martel Calibrated Voltage Source to INPUT 1 on the **BNC to DB9** module. Note: use negative polarity.
- Record 15 seconds of data for each input voltage.

Right two from front.

- - 4 V = -3.997..... volts and GPS Start Time 1252 167 180
 - - 2 V = -1.997..... volts and GPS Start Time 1252 167 262
 - 0 V = -0.0007..... volts and GPS Start Time 1252 167 289

Record Optical Follower Servo settings

- Offset: 4.00..... volt
- Gain: 41.00..... dB
- OFS PD: -3.933..... volt

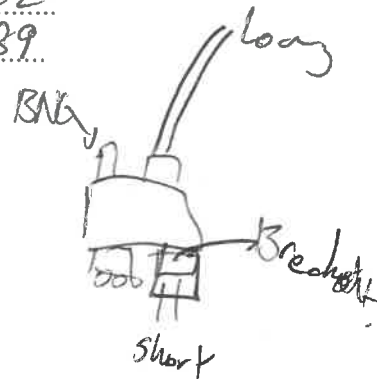

Record Working Standard temperature

- Measure sensor voltage at Interface Module WS PD MON BNC output using a voltmeter.

- GPS time: 1252 167 529
 - WS PD MON: 2.586..... volts

- Break 1k6
Clips from 1k6 to
BNC.

- Convert to degrees Celsius by dividing by 0.90909, multiplying by 100, and subtracting 273.15.

- WS PD temperature:deg. Celsius.
 - WS PD temperature:deg. Fahrenheit.

Time Series Measurements

- Connect Pcal Blue Box PD MON output to INPUT 1 on the **BNC to DB9** module.

- wrap cable around pole.
 Measurement 1:

- Block the OUTER beam with a razor blade dump in the Tx module.
- Place the WS in the INNER beam in the Tx module.

LIGO-T1500062-v8

WS in the INNER beam in the Tx module.			
GPS Times		Readings from MEDM screen	
Start Time #1	1252 167 661	TxPD	2.713 volt
Duration	240 seconds	WSPD	-2.423 volt
End Time #1	1252 167 901	OFSPD	-3.933 volt

Measurement 2:

- Move the block to the INNER beam in the Tx module.
- Move the WS to the OUTER beam in the Tx module.

WS in the OUTER beam in the Tx module.			
GPS Times		Readings from MEDM screen	
Start Time #2	1252 168 104	TxPD	2.714 V
Duration	240 seconds	WSPD	-2.417 V
End Time #2	344	OFSPD	-3.933 V

Measurement 3:

- Leave the WS in the OUTER beam in the Tx module with the INNER beam blocked.
- Close the shutter in the Tx module.

WS in the OUTER beam in the Tx module. Shutter CLOSED.			
GPS Times		Readings from MEDM screen	
Start Time #3	1252 168 364	TxPD	0.0058
Duration	60 seconds	WSPD	-0.0002
End Time #3	424	OFSPD	-0.0094

Measurement 4:

- Leave the block in the INNER beam in the Tx module.
- Open the shutter in the Tx module.
- Replace the Rx sensor with the WS in the Rx module.

WS in the Rx module. INNER beam blocked in the Tx module.			
GPS Times		Readings from MEDM screen	
Start Time #4	1252 168 859	TxPD	2.713
Duration	240 seconds	WSPD	-2.386
End Time #4	9 099	OFSPD	-3.933

Measurement 5:

- Move the block to the OUTER beam in the Tx module.

WS in the Rx module. OUTER beam blocked in the Tx module.			
GPS Times		Readings from MEDM screen	
Start Time #5	1252 169 148	TxPD	2.714
Duration	240 seconds	WSPD	-2.392
End Time #5	388	OFSPD	-3.932

subsubsection*Measurement 6:

- CLOSE the shutter in the Tx module.

WS in the Rx module. Shutter CLOSED in the Tx module.			
GPS Times		Readings from MEDM screen	
Start Time #6	1252 169 452	TxPD	0.0056
Duration	60 seconds	WSPD	-0.0004
End Time #6	512	OFSPD	-0.009

Measurement 7:

- OPEN the shutter in the Tx module.
- Leave the beam block in the OUTER beam in the Tx module.
- Replace the WS with the Rx sensor in the Rx module.

OUTER beam blocked in the Tx module. Rx sensor in the Rx module.			
GPS Times		Readings from MEDM screen	
Start Time #7	1252 169 800	TxPD	2.715
Duration	240 seconds	RxPD	1.724
End Time #7	7 0040	OFSPD	-3.933

Measurement 8:

- Move the beam block to the INNER beam in the Tx module.

INNER beam blocked in the Tx module. Rx sensor in the Rx module.			
GPS Times		Readings from MEDM screen	
Start Time #8	1252 170 075	TxPD	2.714
Duration	240 seconds	RxPD	1.717
End Time #8	315	OFSPD	-3.932

Measurement 9:

- CLOSE the shutter in the Tx module.

Shutter CLOSED in the Tx module. Inner beam blocked.			
GPS Times		Readings from MEDM screen	
Start Time #9	1252 170 336	TxPD	0.0057
Duration	60 seconds	RxPD	0.0009
End Time #9	96	OFSPD	-0.0101

When measurements are finished.

- Remove the beam block from the INNER beam in the Tx module.
- OPEN shutter in the Tx module.
- Set shutter control to *Remote* on interface module.
- Re-enable excitations on the Pcal MEDM screen (if applicable).
- Check that ETM is pointed properly. Check Pcal beam locations at the Rx sensor (photograph spot locations on white card).

LIGO-T1500062-v8

- Analyze the data and upload results to the SVN.
- Make aLog entry, append image of beam spots at the Rx module, add pointer to measurements results in the SVN.

Data Retrieval and Ratio and Force Coefficient calculations

Data Acquisition, Plots and Report

1. Make sure the SVN directory
`/ligo/svncommon/CalSVN/aligocalibration/trunk/Projects/PhotonCalibrator/scripts/pcalEndstation`
is installed on the machine where you are working.
Change to this directory.
Note: A description of the Photon Calibrator SVN directory structure and instructions for checking out the whole Pcal repository can be found in LIGO-T1500095.
2. Execute the command `svn update` to make sure you have the latest versions of the scripts.
3. Execute the command `kinit albert.einstein@LIGO.ORG` to establish connection to the external data server. This will be required for obtaining data later.
4. Navigate to `/ligo/svncommon/CalSVN/aligocalibration/trunk/Projects/PhotonCalibrator/scripts/pcalEndstation`
5. Open the `endstationPcalAnalysis.m` matlab script.
 - (a) Enter the appropriate ifo. location and arm and date code.
 - (b) Make sure the `workingcopy_location` in the Matlab file `pcalworkingcopylocation.m` has the appropriate path to the PhotonCalibrator directory, e.g. `/ligo/svncommon/CalSVN/aligocalibration/trunk/Projects/`
 - (c) If you are running with this calibration data for the first time, make sure that the `rerun` parameter is set to 0.
 - (d) Enter the GPS times of the various measurements between lines 40 and 60.
 - (e) run `endstationPcalAnalysis.m`
6. It will create a folder named `tDYYYYMMDD` in the appropriate end-station-specific directory, e.g. for LHOY
`/ligo/svncommon/CalSVN/aligocalibration/trunk/Projects/PhotonCalibrator/measurements/LHO_EndY`
7. This will also fetch the data from the server and write them into `.txt` files in the `DYYYYMMDD` directory.
8. It also generates plots of the times series of the ratios and saves the plots to the same directory, in addition to displaying them on the screen. Make sure the plots are satisfactory before closing them.
9. It will generate the files, `DYYYYMMDD_Ratio.mat` and `DYYYYMMDD_Results.mat`, that contain the calibration results, and save them in the same directory.

10. Finally it publishes a report LHOY_PD_trend.pdf in the output tDYYYYMMDD directory, and copies over the scripts used for analysis to the sub-directory tDYYYYMMDD/scripts
11. Use the times series plots and the trend document to assess the quality of the recently acquired results.
12. If the data appears to be valid, change the name of the directory from tDYYYYMMDD to DYYYYMMDD. If the data is corrupted in some way (e.g. a large glitch in the data or a data dropout), change the directory name to xDYYYYMMDD. This vetoes the directory from trend documents published at a later date.

Force Coefficient Calculation

1. If the data was not vetoed, then navigate to the directory containing the scripts for making force coefficient trends, in /ligo/svncommon/CalSVN/aligocalibration/trunk/Projects/PhotonCalibrator/pcalCalFactor/scripts
2. Open pcalFCPublish.m, and change to the appropriate date_code. Run the script by passing the appropriate ifo-arm arguments For example for LHOY, run pcalFCPublish('LHOY, '').
3. It creates a force coefficient trend document back in the measurements directory. For LHOY, it will be named pcalFCReport_LHOY.pdf.

Commit the new files to the SVN

1. In the terminal window, execute the following commands (Refer to T1500095 for details) from the
.../PhotonCalibrator/measurements/end_station
folder (Refer to T1500095 for details.)
 - svn add DYYYYMMDD
 - svn commit -m "message: including comment on veracity of data and analysis results by person committing the directory to the SVN"
 - svn update